



3800 N. Mill Road Vineland, NJ 08360 Phone: 800.843.1794

Fax: 800.922.4361

AF-0300 Vacuum Cart Systems

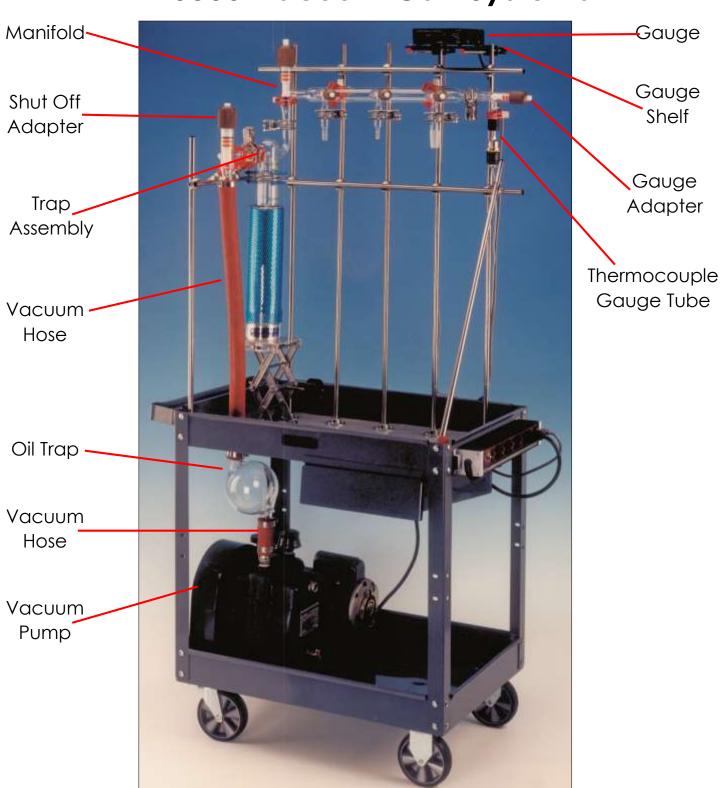

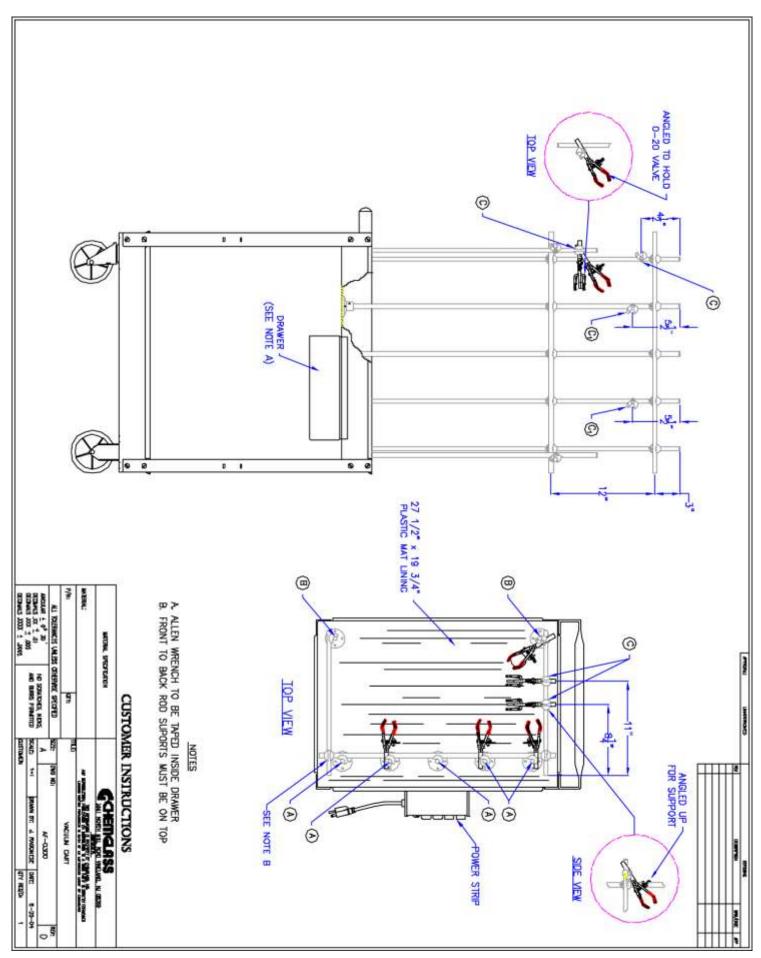

Assembly Instructions for AF-0300 Vacuum Cart Systems

Your cart contains the following stainless steel rods:

5 each: 39"OAL2 each: 26.5"OAL2 each: 29"OAL2 each: 16.25"OAL

To Assemble:

- Insert 5 each 39"rods in support feet (A)
 (DO NOT TIGHTEN ALLEN SCREWS INTO SUPPORT FEET NOW)
- Insert 2 each 29"rods in support feet (B)
 (DO NOT TIGHTEN ALLEN SCREWS INTO SUPPORT FEET NOW)
- 3. Install 2 each 26.5" rods horizontally on the uprights (as per blueprint)
- 4. Install 2 each 16.25" rods horizontally. Secure between front and back supports. (See note 1 on blueprint)
- 5. Tighten all Allen screws on the support feet.

<u>Mounting of Glassware on Lattice System:</u>

The enclosed blueprint drawing shows the approximate dimensions for mounting the glassware on the lattice using the clamps. The bosshead locations are marked as "C".

To ensure that the manifold and trap assemblies are square and level, the back of the clamp rods are to be even with each other, marked as C^1 .

The rubber vacuum tubing supplied with the cart fits snugly within the black rubber grommet. When installing the tubing, the grommet must be removed from the cart and slid over the tubing (above cart). After the tubing is tightened to the main shut off and oil trap (or pump), the grommet can be reinstalled into the opening around the tubing.

If you have any questions, please contact us:



3800 North Mill Road, Vineland, NJ 08360 US & Canada Phone: (800) 843-1794 International: (856) 696-0014 US & Canada Fax: (800) 922-4361 International: (856) 696-9102

^{*}Allen wrench is in the drawer of the cart.