


CG-1950 & CG-1951 Benchtop Support Stands

Single, Double, & Mobile

Each Kit Contains The Following:

#	Qty	Description
A	1	Reactor Support Frame Base
B	2	S.S. Pipe for Uprights
C	2	Crossbar for Motor/Reactor Support
D	1	Single Reactor Base Plate
E	1	Dual Reactor Base Plate
F	2	Reactor Support Rod for 2L to 5L
G	2	Reactor Support Rod for 250mL to 1L with Velcro
H	1	120mm ID Cork Ring fits 2L to 5L
I	2	90mm ID Cork Ring fits 300mL to 1L
J	2	60mm ID Cork Ring fits 250mL

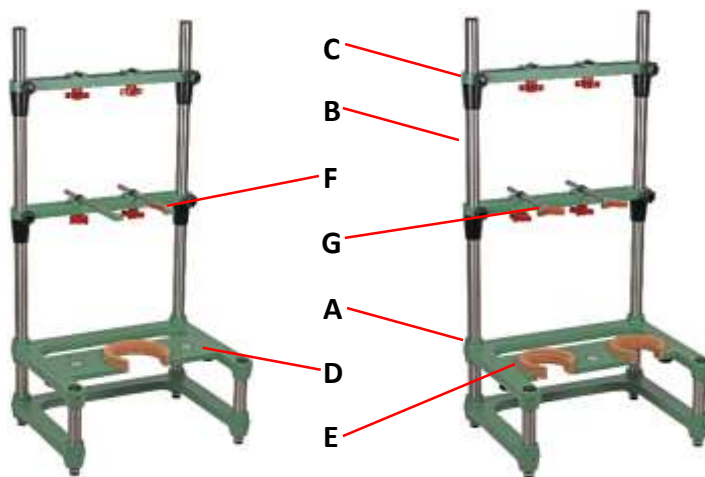

Assembly Instructions:

1. Remove the support plates from the frame base.
2. Turn the base over on its side. Install the uprights by sliding the S.S. pipe through the bottom of the fittings in the base. Tighten the Allen screws. Note: Leveling feet should be at the bottom.
3. With the stand in the upright position- slide both crossbars over the uprights and position as shown on the diagram. The red knobs should be on the bottom.
4. Tighten the black knobs on each end to hold them in place. The positions will have to be adjusted depending on the glassware used in the stand.
5. Install the support rod(s), cork ring(s) and the base plate that corresponds with the reactor size that you are using.

