


CONTACT YOUR IT SPECIALIST FOR ALL THE INFORMATION ABOUT YOUR WIFI CONNECTION

1. ACTIVATE ERMES ACCOUNT

- **Create** Your VELP Account or **login** if you already have it.
- Access to your profile page and select “**Configure Your VELP Ermes Account**”. Accept Terms and Conditions and **Activate your Ermes account**.
- An email will be send to the email address that you have insert during the registration containing your **ID, Password** and a **link** to proceed to VELP Ermes authentication.
- At the first login on VELP Ermes a new password will be request: *remember to change it!*

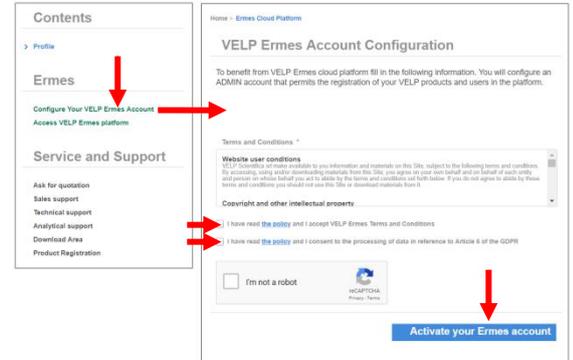


2. CONNECT TO YOUR WIFI CONNECTION

- Turn on your AREX-6 Connect PRO, select the VELP_AREX6_(serial number) available on the Wi-Fi list, in order to connect the instrument.

Note If the instrument doesn't appears in the list or it was already configured with another WiFi, follow the **RESET procedure**:

1. Push for 3 seconds both knobs
2. Push the central button until is displayed STAR AP
3. Push and Rotate the right knob to select YES
4. Push the right knob to confirm
5. Select VELP_AREX6_(serial number) available on the Wi-Fi list, in order to connect the instrument.

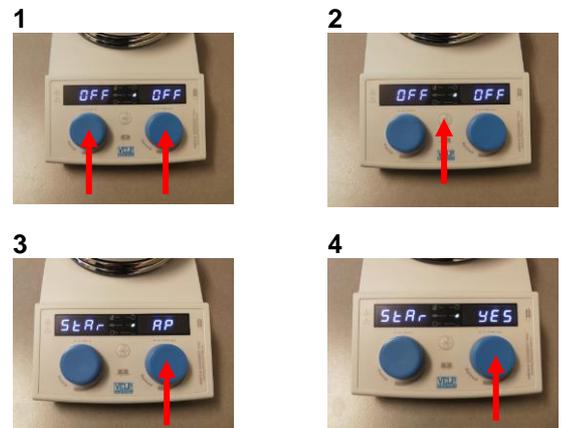

- Open a browser from your device, insert the address **192.168.0.50** and set the parameters required to connect the instrument to your Wi-Fi and press on Push to confirm. **As Passkey, insert: velp**



EXAMPLE:


PassKey: velp
SSID: (your WiFi name)*
Pwd: (your WiFi password)*
Auth: Wpa & WPA2 Personal

Press on **Push** to confirm and wait until the “This site can't be reached” is displayed on your screen.

*If necessary, please contact your IT administrator.

- Wait until the Wi-Fi LED Icon stop blinking.

NOTE: If the LED is still blinking after 2 minutes, please try the procedure again and check if connection parameters are correct.

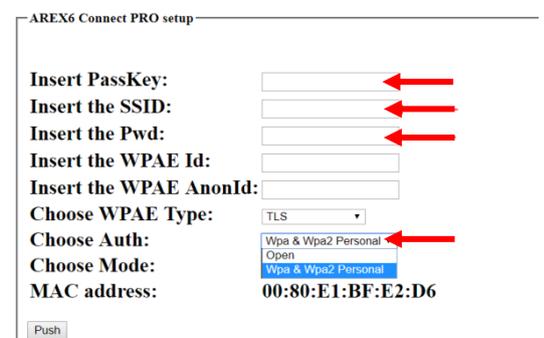

3. REGISTRATION ON ERMES PLATFORM

1

- Push for 3 seconds both knobs
- Push the central button until is displayed OTC
- Push the right knob in order to see - - - - on display




2

Login on VELP Ermes with your ID and Password received by email, and select "**Device Registration**".



3


Insert the AREX-6 Connect PRO **serial number**, **product code** (you can find these information on the silver label on the back of it) and **purchase date**. Press Next.

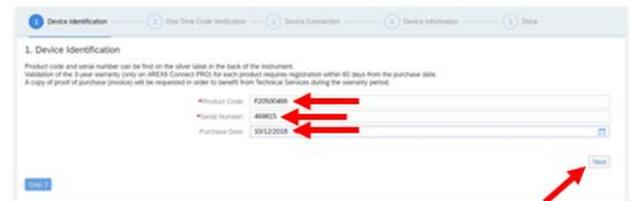

4


On the display of AREX-6 Connect PRO will appears a One Time Code (OTC) of four digits. This number must be insert for security reasons. Press Next and wait for connection.

NOTE: If the OTC doesn't appear on AREX-6 Connect PRO, please contact your IT administrator and perform again the step **Connect to your WiFi connection**

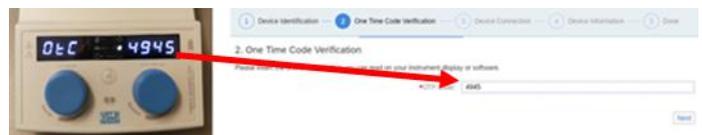


5

Give a **name** to your AREX-6 Connect PRO and inset the **address** where your instrument is located. Press Next

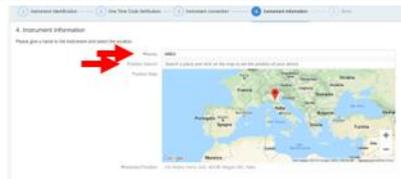


6

From the homepage select "**Laboratory Management**", then add the AREX-6 Connect PRO to an existing laboratory by pressing edit and "+", or create a new one and add the AREX-6 Connect PRO by pressing "+", then add users to the laboratory.

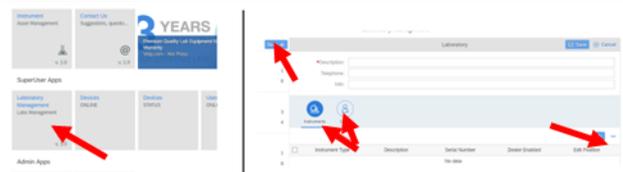

7

Now from "**Instruments**" it is possible to monitor and control your connected AREX-6 Connect PRO.

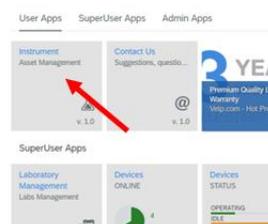

WELCOME TO THE SMART LAB!

Elevating laboratories to the cloud