**Instruction Manual**  
**Manuale di istruzioni**  
**Manuel d'instructions**  
**Manual de instrucciones**  
**Bedienungsanleitung**  
**指导手册**

**OHS 200 DIGITAL Overhead stirrers**

F20100490, F20110490

**OHS 100 DIGITAL Overhead stirrers**

F20100491, F20110491

**OHS 60 DIGITAL Overhead stirrers**

F20100492, F20110492

**OHS 40 DIGITAL Overhead stirrers**

F20100493, F20110493

**OHS 20 DIGITAL Overhead stirrers**

F20100494, F20110494

**General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息**

Before using the unit, please read the following instruction manual carefully.  
Prima dell'utilizzo dello strumento si raccomanda di leggere attentamente il seguente manuale operativo.  
Avant d'utiliser l'instrument, il est recommandé de lire attentivement le présent manuel d'instructions.  
Antes de utilizar el instrumento, le recomendamos que lea con atención el siguiente manual de funcionamiento.  
Bitte lesen Sie vor Inbetriebnahme des Geräts diese Bedienungsanleitung sorgfältig durch  
在使用本装置之前，请仔细阅读以下使用说明书。

Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.  
Non smaltire l'apparecchiatura come rifiuto urbano, secondo quanto previsto dalla Direttiva 2002/96/CE.  
Ne pas recycler l'appareil comme déchet solide urbain, conformément à la Directive 2002/96/CE.  
No tirar el aparato en los desechos urbanos, como exige la Directiva 2002/96/CE.  
Dieses Gerät unterliegt der Richtlinie 2002/96/EG und darf nicht mit dem normalen Hausmüll entsorgt werden.  
根据 EEC 指令 2002/96/CE，请不要将本设备作为城市垃圾处理。

**This unit must be used for laboratory applications indoor only.** The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with not specified accessories, product's safety may be compromised.

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**This unit has been designed and manufactured in compliance with the following standards:**

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Electrical equipment for laboratory use

**UL 61010-1**

General requirement - Canadian electrical code

**CAN/CSA-C22.2 No.61010-1**

VELP reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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The values indicated on the rating plate of the instrument must correspond to those of the power supply.

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Position the instrument on a flat surface, with a distance from the wall of 30 cm (at least).

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Safe working conditions are ensured only when the accessories described in the dedicated chapter are used.

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Do not use with explosive or dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie or to stir harmful liquids prior using protective measures according to the safety standards of the processed products and/or in force in the laboratories including personal protective equipment and the presence of an extraction hood which ensures at least 10-fold air change in accordance with the standards EN 14175 and DIN 12924.

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It is dangerous to run the unit with the stirring blade turning in free air. Always place the stirring shaft in the receptacle before turning the unit on.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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The solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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Switch off the the stirring before opening chuck locking ring. Switch off the instrument before removing chuck.

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OHS overhead stirrers with electronic speed control, brushless motor, and advanced safety features are able to satisfy the most difficult laboratory applications in terms of viscosity and volume. The new chuck ensures higher safety for the operator and allows to use a passing rod of up to 8.5mm in diameter.

Figure 2. OHS 200 Digital

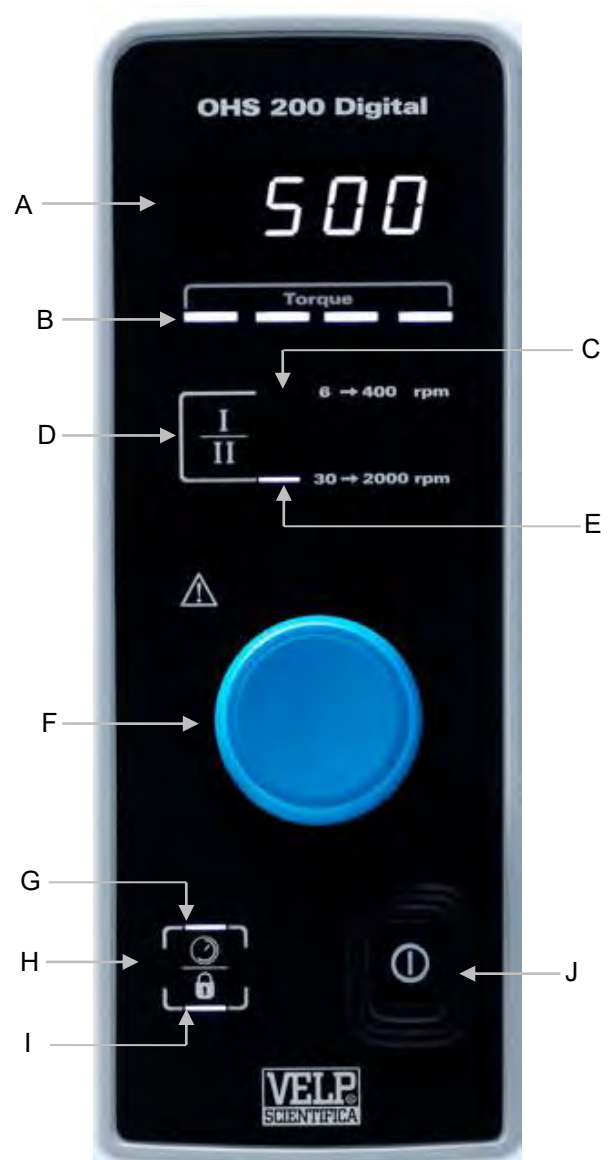
Figure 1. OHS 200 Digital - Frontal view

- A 7 Segment Display LED
- B Torque indicator bar
- C Low speed LED (only OHS 200)
- D Gear key (only OHS 200) – Timer key for other models
- E High speed LED (only OHS 200) – Timer LED for other models

- F Speed control knob
- G Timer LED (for OHS 200) - Upper lock LED for other models
- H Timer key (for OHS 200) – Lock key for all models
- I Lower Lock LED
- J Main switch
- K Smart chuck

## 2. Assembly and installation

- Unpacking
  - Check the integrity of the unit after unpacking.
- The box includes
  - OHS (XXX) digital overhead stirrer
  - Power supply cable
  - Instruction manual
  - Allen key and allen head screw
  - Shaft support overhead stirrer
- Installation
  - Place the unit on non-flammable surface
  - Fix the shaft support to the instrument with the allen head screw
  - Fasten the unit to the support rod (A00000366 or A00000369) using the double clamp (A00001301)
  - Secure the receptacle using the ribbon clamp (A00001302)
  - Slide the stirring shaft into the chuck and tight it
  - Make sure that the rating value of the instrument corresponds to the one of the power supply
  - Ensure that the socket provided, with grounding, is compliant with the current safety norms and easy to reach.

## 3. Working

<b>Commissioning</b>	<ul style="list-style-type: none"> <li>➤ Switch on the instrument through the main switch (J)</li> <li>➤ Display (A) shows software version, last set point value and OFF (if "Mode" is set on Stop, see chapter 0).</li> </ul>
<b>Stirring</b>	<ul style="list-style-type: none"> <li>➤ Set the operating speed range by clicking Gear key (D) (Only OHS 200).</li> <li>➤ Adjust speed set point by turning the speed control knob (F).</li> <li>➤ Click the knob to start stirring.</li> <li>➤ Speed increases until set point achievement.</li> <li>➤ A microprocessor ensures constant speed even when the viscosity changes (counter-reaction).</li> <li>➤ Switch off the stirring by clicking the knob.</li> </ul>
<b>Timer</b>	<ul style="list-style-type: none"> <li>➤ Click Timer key (D or H). HH:MM appears on display.</li> <li>➤ Rotate speed control knob (F) to set the desired value. Click to confirm.</li> <li>➤ Once timer is set, timer LED lights up:               <ul style="list-style-type: none"> <li>- (G) for OHS 200</li> <li>- (E) for other models</li> </ul> </li> <li>➤ The LED lights continuously until the instrument starts to stir. When the instrument stirs, the LED flashes.</li> <li>➤ It's possible to visualize (or modify) countdown in each moment during the analysis by clicking timer key.</li> <li>➤ When countdown finishes, "End" is visualized on the display.</li> </ul>
<b>Gear (only for OHS 200)</b>	<ul style="list-style-type: none"> <li>➤ Set the operating speed range by clicking Gear button (D).</li> <li>➤ The gear set has its relative LED (C or E) lighted on.</li> <li>➤ It's possible to select between two gears:               <ul style="list-style-type: none"> <li>- <b>Gear I:</b> Low speed (<b>6 – 400 rpm</b>) High torque (<b>till 200 Ncm</b>).</li> <li>- <b>Gear II:</b> High speed (<b>30 – 2000 rpm</b>) Low torque (<b>till 40 Ncm</b>).</li> </ul> </li> </ul>
<b>Torque</b>	<ul style="list-style-type: none"> <li>➤ The torque indicator bar (B), made by four LEDs, shows the intensity of the torque applied</li> </ul>
<b>Lock</b>	<ul style="list-style-type: none"> <li>➤ Holding the Lock key (H) for 3 seconds it's possible to lock the control panel.</li> <li>➤ Unlock the control panel by holding key H for 3 seconds.</li> <li>➤ When the instrument is locked,               <ul style="list-style-type: none"> <li>- LED (I) lights on for OHS 200</li> <li>- LED (G) and (I) light on for other models</li> </ul> </li> </ul>

## 4. External Connections

- |            |   |
|------------|---|
| <b>USB</b> | <ul style="list-style-type: none"> <li>➤ Each model has USB connection for software upgrading by PC. Software version, dedicated program to be installed on the PC and software installation guide must be requested by e-mail to <a href="mailto:service@velp.it">service@velp.it</a></li> </ul> |
|------------|---|

## 5. Menu

Hold the control knob (F) for 3 seconds to enter into the setting mode when stirring is off.

Turning the control knob (F) it is possible to pass from one of the following parameters to the next one. Click the knob to modify:

Parameter shown on display	Default value		Range	Description
NOdE	StOP		StOP - run	Set starting mode: <ul style="list-style-type: none"> <li>➤ Stop: when the instrument is switched on through the main switch, OFF is on display</li> <li>➤ Run: when the instrument is switched on, it restarts to work with the last set points set</li> </ul>
LINT	OHS 200	400 (I) 2000 (II)	100 – 400 100 - 2000	Set the maximum value of speed set point speed
	OHS 100	1300	100 – 1300	
	OHS 60	2000	100 – 2000	
	OHS 40	2000	100 - 2000	
	OHS 20	2000	100 - 2000	
COun				Running time in H. If “.” Is showed before the number, the running time is in Day.
UPGr				To install a new software version
rSEt	nO		YES - nO	To restore default setting parameters
CAI	nO		YES - nO	To reset torque zero value

To leave settings mode wait 10 seconds without touching any knob or click the lock key.

## 6. SmartChuck™

Smart chuck (K) allows to fix stirring shaft with one hand.

- Open the chuck: From work position (Figure 3), turn Locking Ring 90° to the left (Figure 4). Pull Locking Ring down (Figure 5).
- Close the chuck: Push Locking Ring up (Figure 6) and turn it 90° to the right.

Figure 3. Working position

Figure 4. Intermediate position

Figure 5. Open position

Figure 6. Intermediate position

## 7. Error messages

When the display shows an error message, the stirring function stops automatically.

Error code	Cause
AL1	Motor doesn't start stirring
AL2	High internal motor temperature
AL3	Motor overload
AL4	High driver temperature
AL5	Safety relay fault

To remove the error message, disconnect the instrument from the power supply.

If alarm persists on the display, please contact VELP Scientifica's technical service department. [service@velp.it](mailto:service@velp.it)

## 8. Maintenance

<b>Maintenance</b>	<ul style="list-style-type: none"> <li>➤ No routine or extraordinary maintenance is necessary.</li> <li>➤ Repairs must be carried out by authorized Velp personnel only.</li> <li>➤ Instrument must be transported in its original packaging any indications present on the original packaging must be followed (e.g. palletized).</li> <li>➤ It is the responsibility of the user, to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.</li> </ul>
<b>Cleaning</b>	<ul style="list-style-type: none"> <li>➤ Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent.</li> </ul>

## 9. Technical data

Models		F20100490, F20100491, F20100492, F20100493	F20110490, F20110491, F20110492, F20110493	
Power supply		230 V / 50-60 Hz (+/-10%)	115V / 60 Hz (+/-10%)	
Dimensions (WxHxD)	OHS200	90x315x235 mm (3.54x12.40x9.25 in)	90x315x235 mm (3.54x12.40x9.25 in)	
	OHS 100 – OHS60 – OHS40 – OHS20	90x285x235 mm (3.54x11.22x9.25 in)	90x285x235 mm (3.54x11.22x9.25 in)	
Weight	OHS200	4,6 kg (10,14 lb)	4,6 kg (10,14 lb)	
	OHS100	4,1 kg (9,04 lb)	4,1 kg (9,04 lb)	
	OHS60 – OHS40 – OHS20	4,1 kg (9,04 lb)	4,1 kg (9,04 lb)	
Power input		190 W	190 W	
Construction material (structure)		Aluminum	Aluminum	
Working in continuous		Admitted	Admitted	
Settable restart modality		Stop or work	Stop or work	
Noisiness		<< 60 dBa	<< 60 dBa	
Environmental temperature admitted		+5...+40 °C	+5...+40 °C	
Storage temperature admitted		-10...+60 °C	-10...+60 °C	
Max humidity		80%	80%	
Level of electrical protection CEI EN60529		IP 54	IP 54	
Overvoltage category		II	II	
Pollution degree CEI EN61010-1		2	2	
Max altitude		2000 m	2000 m	
<b>Stir</b>	Stirring capacity	OHS200	100 l H <sub>2</sub> O	100 l H <sub>2</sub> O
		OHS100	100 l H <sub>2</sub> O	100 l H <sub>2</sub> O
		OHS60	40 l H <sub>2</sub> O	40 l H <sub>2</sub> O
		OHS40	25 l H <sub>2</sub> O	25 l H <sub>2</sub> O
		OHS20	25 l H <sub>2</sub> O	25 l H <sub>2</sub> O
	Programmable speed range	OHS200	6-400rpm (I) – 30-2000rpm (II)	6-400rpm (I) – 30-2000rpm (II)
		OHS100	30-1300rpm	30-1300rpm
		OHS60	30-2000rpm	30-2000rpm
		OHS40	30-2000rpm	30-2000rpm
		OHS20	30-2000rpm	30-2000rpm
Motor type		BLDC	BLDC	
Speed selection		1 rpm step	1 rpm step	
Stirring alarm		Motor fault	Motor fault	
Motor rating output		150 W	150 W	
<b>Torque</b>	Max torque admitted	OHS200	200 Ncm (I) – 40 Ncm (II)	200 Ncm (I) – 40 Ncm (II)
		OHS100	100 Ncm	100 Ncm
		OHS60	60 Ncm	60 Ncm
		OHS40	40 Ncm	40 Ncm
		OHS20	20 Ncm	20 Ncm
<b>Counters</b>	Motor counter	Working hours	Working hours	

## 10. Accessories / Spare parts

A00000366	Support rod and base OHS	A00001307	Stirring shaft with propeller
A00000369	Telescopic support rod and base OHS	A00001308	Stirring shaft with paddle, 6 holes
A00000391	OHS/ControllerSoft	A00001309	Stirring shaft with turbine
A00001300	Support rod and base	A00001310	Stirring shaft with turbo propeller
A00001301	Double clamp	A00001311	Stirring shaft with anchor
A00001302	Ribbon clamp	A00001312	Chuck guard OHS
A00001304	Stirring shaft with floating blade		
A00001305	Stirring shaft with folding blade	10007175	Knob 35D blu
A00001306	Stirring shaft with fixed blade		

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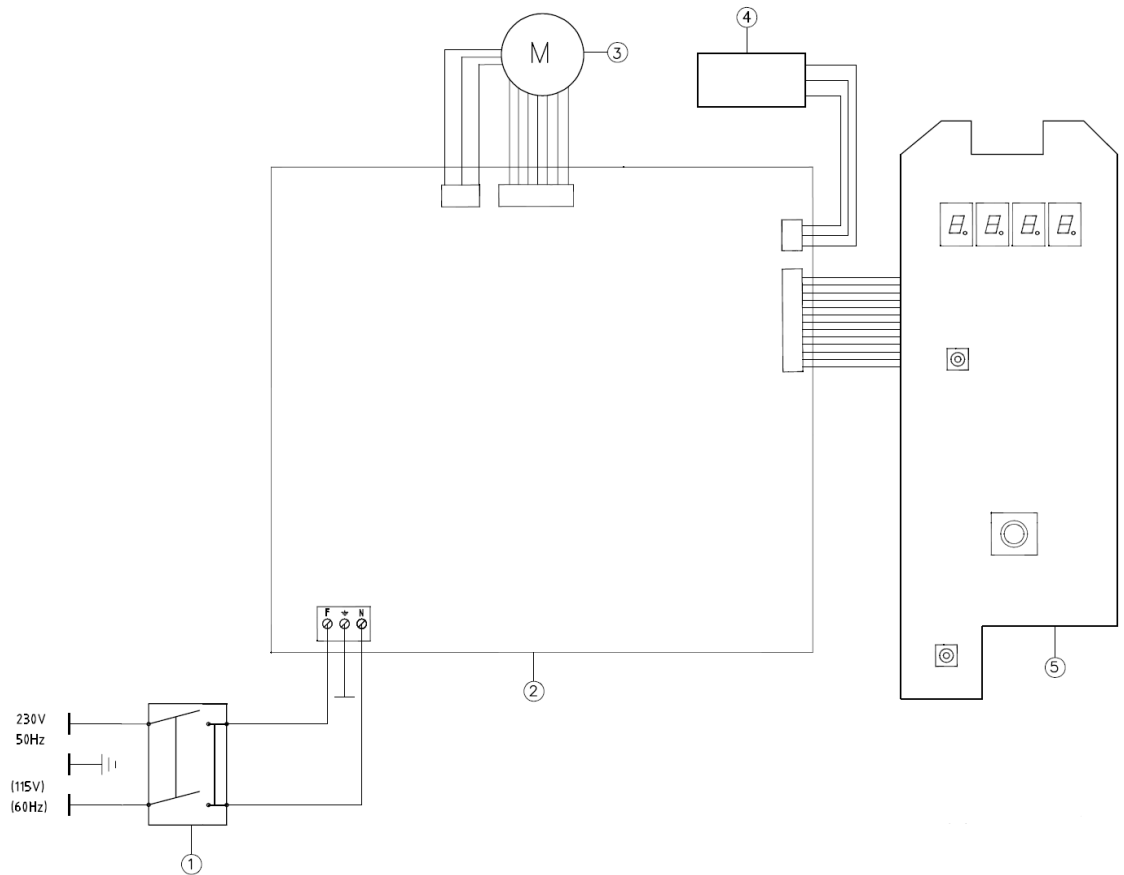
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1	Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use - Part 1: General Requirements
EN 61010-2-051	Safety requirements for electrical equipment for measurement, control and laboratory use - Part 2-051: Particular requirements for laboratory equipment for mixing and stirring
EN 61326-1	Electrical equipment for measurement, control and laboratory use - EMC requirements -
2015/863/EU (RoHS III)	Restriction of the use of certain hazardous substances
2012/19/EU (WEEE)	On waste electrical and electronic equipment

and satisfies the essential requirements of the following directives: / e soddisfa i requisiti essenziali delle direttive: / et qu'il satisfait les exigences essentielles des directives: / y cumple con los requisitos esenciales de las directivas: / und den Anforderungen folgender Richtlinien entspricht: / 并满足以下指令的基本要求。

- Machinery directive 2006/42/EC / Macchine 2006/42/CE / Machines 2006/42/CE / Máquinas 2006/42/CE / Maschinen 2006/42/EG / 机械指令 2006/42/EC
- Electromagnetic compatibility directive 2014/30/EU / Compatibilità elettromagnetica 2014/30/EU / Compatibilité électromagnétique 2014/30/EU / Compatibilidad electromagnética 2014/30/EU / Elektromagnetische Verträglichkeit 2014/30/EU / 电磁兼容性指令 2014/30/EU
- plus modifications / più modifiche / plus modifications / más sucesivas modificaciones / in der jeweils gültigen Fassung / 加上变化

## 1. Declaration of conformity <sup>UK</sup><sub>CA</sub>

We, the manufacturer VELD Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use  
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -  
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment  
Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

## Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

### Analytical instruments

Elemental Analyzers  
Digestion Units  
Distillation Units  
Solvent Extractors  
Fiber Analyzers  
Dietary Fiber Analyzers  
Oxidation Stability Reactor  
Consumables

### Laboratory Equipment

Magnetic Stirrers  
Heating Magnetic Stirrers  
Heating Plates  
Overhead stirrers  
Vortex mixers  
Homogenizers  
COD Thermoreactors  
BOD and Respirometers  
Cooled Incubators  
Flocculators  
Overhead Shakers  
Turbidimeter  
Radiation Detector  
Open Circulating Baths  
Pumps

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VELP Scientifica Srl