Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
指导手册

OV 625 Digital Dispenser

F20900470, F20910470

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 一般信息

Before using the unit, please read the following instruction manual carefully.
Prima dell'utilizzo dello strumento, si raccomanda di leggere attentamente il seguente manuale operativo.
Avant d'utiliser l'instrument, il est recommandé de lire attentivement le présent manuel d'instructions.
Antes de utilizar el instrumento, le recomendamos que lea con atención el siguiente manual de funcionamiento.
Bitte lesen Sie vor Inbetriebnahme des Geräts diese Bedienungsanleitung sorgfältig durch
在使用本装置之前，请仔细阅读以下使用说明书。

Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.
Non smaltire l'apparecchiatura come rifiuto urbano, secondo quanto previsto dalla Direttiva 2002/96/CE.
Ne pas recycler l'appareil comme déchet solide urbain, conformément à la Directive 2002/96/CE.
No tirar el aparato en los desechos urbanos, como exige la Directiva 2002/96/CE.
Dieses Gerät unterliegt der Richtlinie 2002/96/EG und darf nicht mit dem normalen Hausmüll entsorgt werden.
根据 EEC 指令 2002/96/CE，请不要将本设备作为城市垃圾处理。

This unit must be used for indoor laboratory applications only. The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a way not specified by the manufacturer or with non-specified accessories, the product's safety may be compromised.

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This unit has been designed and manufactured in compliance with the following standards:

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Electrical equipment for laboratory use
General requirement - Canadian electrical code

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safe substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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The solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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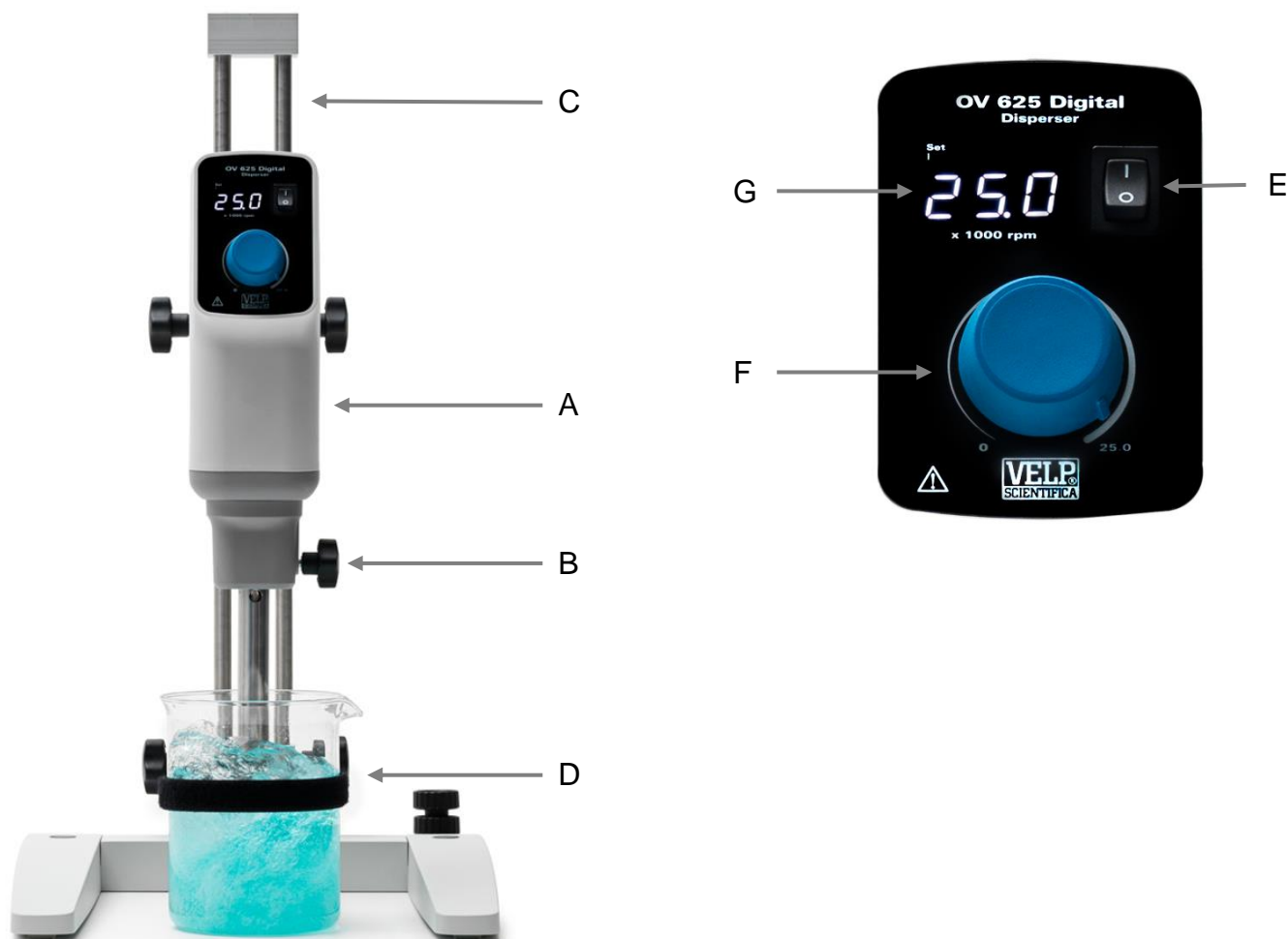
The OV 625 Digital Disperser, equipped with electronic speed control system and brushed motor, it is the ideal solution for dispersing, homogenizing, mixing and crushing samples of biological materials (cells, animal and plant tissues), pharmaceutical, cosmetic and food products.

The term dispersion means the subdivision and distribution of a solid or fluid in a liquid phase that is not completely miscible with the other phase.

The dispersing tool is mainly composed of the rotor and the stator. The high number of revolutions of the rotor allow the tool to axially suck the sample into the dispersion head and then push it radially through the slots between the rotor and the stator. Normally, the processing time is a few minutes to obtain the desired fineness, longer processing times only slightly improve the fineness of the sample but significantly increase its temperature.

	Description	Function
A	OV 625 Digital	D20 Drive unit
B	Dispersing tool fixing knob	To fix the dispersing tool into disperser
C	Mini H stand	To fix the disperser
D	Fixing clamp	To fix the receptacle with the sample
E	Main switch	To turn the disperser on
F	Control knob	Rotate to start/stop the dispersing process, adjust the speed
G	3-digit LED display	To show the speed value (\div 1000 rpm)

2. Assembly and installation

- Unpacking
 - Check the integrity of the unit after unpacking.
- The box includes
 - OV625 Digital disperser (A)
 - Tool disassembly kit
 - Instruction manual
 - Allen key and allen head screw
 - Support shaft
- Installation
 - Place the Mini H stand (C) (accessory) on a surface as described on safety regulations
 - Fix the support shaft to the instrument with the allen head screw
 - Firmly fix the instrument to the Mini H stand (C)
 - Secure the receptacle with the sample using the fixing clamp (D) supplied to the Mini H stand.
 - Slide the dispersing tool (accessory) into the head of the disperser up to the stop and tight it with the knob (B)
 - Place the dispersing tool into the sample as indicated in Chapter 5
 - Make sure that the rating value of the instrument corresponds to the one of the power supply
 - Ensure that the socket provided, with grounding, is compliant with the current safety norms and easy to reach.

IMPORTANT NOTE

- Check that the support shaft is firmly seated. The vibrations may loosen the allen head screw, for safety reasons, be sure to tighten it periodically.
- Check that the disperser is securely mounted in position before use and at regular periods. The position of disperser can only be changed if the instrument is turned off and disconnected from the power supply

3. Operation

- Power-up
 - Turn the instrument on through the main switch (E)
 - The display (G) shows the software version, the last setpoint value and $\square F F$ (if “ $\square P E$ ” is set on R , see Chapter 0)
- Dispersing
 - Turn the speed control knob (F) to set the speed and start the dispersing process
 - Speed increases until set point achievement
 - A microprocessor ensures constant speed even when the viscosity changes (SpeedServo)
 - Stop the process by turning the control knob to 0 (F)

4. Setting mode

Turn on the instrument using the main switch (E); while the software version is displayed, by turning the speed knob (F) from minimum to maximum at least two times, the hidden parameters are shown.

Rotate the speed knob to scroll through the different menus available.

If the knob is not moved for 2 seconds, the display shows the current value of the selected menu for 5 seconds and then goes back, showing the name of the menu if the knob is not turned. For parameters with the possibility of modification, if the knob is turned (the set LED lights up to indicate that the parameter is being modified), the display shows the parameter values proportionally to the position of the knob.


The new value is saved if you do not turn the speed knob for 4 seconds, the display shows “Eнд” (lighting off the Set led) for 2 seconds before coming back to show the name of the menu.

Parameter	Default value	Range	Description
DPt	R	R - b	Set starting mode: ➤ R: when the instrument is switched on through the main switch, its working state is OFF ➤ b: when the instrument is switched on through the main switch, it restarts to work with the last speed set.
EnE	-	-	Operating time of brushes in hours. If “” is shown in the number, the running time is in days.
EnP	-	-	Operating time of motor in hours. If “” is shown in the number, the running time is in days.

5. Dispersing tools

Insertion of the dispersing tool into the disperser

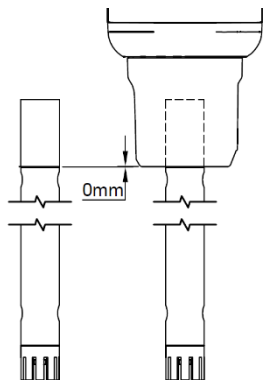Unscrew the knob (B) until its head is no more visible through the hole. Slide the dispersing tool from the bottom into the disperser, up to the stop. The dispersing tool enters with an audible click if you insert it with enough force. It's fitted correctly when the notch on the dispersing tool is aligned with the bottom of the disperser.

Now fix the dispersing tool by tighten the knob (B).

Note: before turning on the unit, make sure that the dispersing tool is inserted in the liquid and that the receptacle is well fixed with the clamp (D).

ABBREVIATIONS

All D20 series tools are allowed (it means with a 20mm diameter head disperser as attachment).

Material: stainless steel AISI 316L.

Head of disperser:

D20 For drive unit with a head of 20mm as diameter

Diameter and type of stator:

S20C Coarse stator diam. 20 mm

S20F Fine stator diam. 20 mm

S25C Coarse stator diam. 25 mm

S25F Fine stator diam. 25 mm

S40C Coarse stator diam. 40 mm (only for mixtures up to 20.000 rpm)

Type of gasket:

P PTFE

Diameter and type of rotor:

R20S Standard rotor diam. 20 mm
R20C Coarse rotor diam. 20 mm
R20E Emulsifying rotor diam. 20 mm
R20M Mixing rotor diam. 20 mm
R25C Coarse rotor diam. 25 mm
R25E Emulsifying rotor diam. 25 mm
R25M Mixing rotor diam. 25 mm

Example of dispersing tool configuration

D20-S20C-P-R20S Velp code: A00000478

AVAILABLE DISPERSING TOOLS

See all dispersing tools (accessories) in Chapter 9 or go to www.velp.com.

Use only the original Velp Scientifica dispersing tools after reading their instruction manual with more information about applications and technical data.

TO OBTAIN BETTER RESULTS

Diameter of the receptacle with the sample

It is better to disperse using a container slightly larger than the dispersing tool in order to obtain a faster and more homogeneous solution.

This will prevent part of the sample from remaining on the walls of the container and therefore not being homogenized.

Keep in mind that the insertion of the tool into the receptacle causes the liquid level to rise.

Position of the dispersing tool

- Place the dispersing tool not in the middle of the receptacle (Figure 1)
- Place the dispersing tool about 1/3 of the sample height from the bottom of the receptacle (Figure 2).

Figure 1

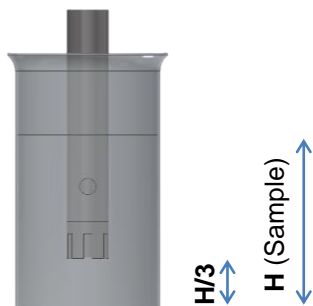
Figure 2

If you use a dispersing tool that is too large for the sample amount, it could suck air.

Dispersing speed

Always start at a low speed, then gradually increase it. This allows to keep control of the process.

Normally the dispersion occurs in 30 – 60 seconds, if this does not happen, check:

- The dispersing tool position in the container
- If the sample is too viscous
- Incorrect dispersing tool for the sample (size, type)

TO NOT REDUCE THE LIFETIME OF THE DISPERSING TOOL AND THE DISPERSER

Never run the dispersing tool without liquid

The dispersing tool requires a liquid sample to lubricate the gasket during machining.

A PTFE gasket is placed between the rotor and stator to prevent the rotor from bending or moving inside the tool. The gasket must be immersed in the liquid sample, otherwise it will burn causing overheating and subsequent damage to the sample and the dispersing tool.

Without the liquid, the PTFE gasket can burn causing damage to the tool and the disperser.

Cleaning and maintenance of dispersing tool

It's necessary to disassemble and wash the dispersing tool after each use, checking the good condition of the gasket and the metal parts and replacing them if necessary. The PTFE gasket is naturally subjected to wear because of the friction with the rotating elements of the tool; therefore it's necessary to periodically replace it in order to guarantee the correct operation of the tool and the dispersing instrument.

Replace the PTFE gasket if the following conditions occur:

- Visible wear of the component (e.g. deformations, thinning),
- Unusual noise and vibration during the operation of the dispersing instrument,
- When the space between the gasket and the internal and external metal parts is excessive.

To disassemble the dispersing tool, use the appropriate tool disassembly kit supplied with the drive unit and read the instruction manual supplied with each tool.

6. Error and warning messages

When the display shows an error message, the motor stops automatically. To restore the functionality of the instrument, proceed as follows:

- Turn the unit off using the main switch
- Use the suggestions described in the column "Solution"
- Turn the unit on

Code	Cause	Solution
RL1	The motor doesn't start or doesn't reach the set speed	<ul style="list-style-type: none"> • Check that there aren't locks in the dispersing tool or structure • Check the operating time of the motor and brushes (chapter 0)
RL2	High internal motor temperature	<ul style="list-style-type: none"> • Check that there aren't obstructions on the air vents for internal ventilation • Check that there aren't locks in the dispersing tool or structure • Reduce speed or load
RL3	Motor overload	<ul style="list-style-type: none"> • Check that there aren't obstructions on the air vents for internal ventilation. • Check that there aren't locks in the dispersing tool or structure • Reduce speed or load
RL5	Problem on the main board	<ul style="list-style-type: none"> • Contact VELP Scientifica's technical service department
RL8	Speed variation too high	<ul style="list-style-type: none"> • Check that there are no locks in the dispersing tool or structure • Check the operating time of the motor and brushes (chapter 0)
RL9	Blackout or plug inserted with the main switch in position I	<ul style="list-style-type: none"> • Enter the plug when the main switch is in position 0

If an error persists on the display, please contact VELP Scientifica's technical service department.

7. Maintenance and cleaning

No routine or extraordinary maintenance is necessary for the unit. The maintenance is foreseen only on the dispersing tools as described in Chapter 5.

Cleaning

Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent. During cleaning operations, prevent water and humidity residues from entering the unit.

Repair

Repairs must be carried out by authorized Velp personnel only.

The transport of the instrument by freight forwarders, couriers or others must be carried out using the original shockproof packaging. Follow any instructions on the original packaging (e.g., palletizing).

It is the responsibility of the user to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

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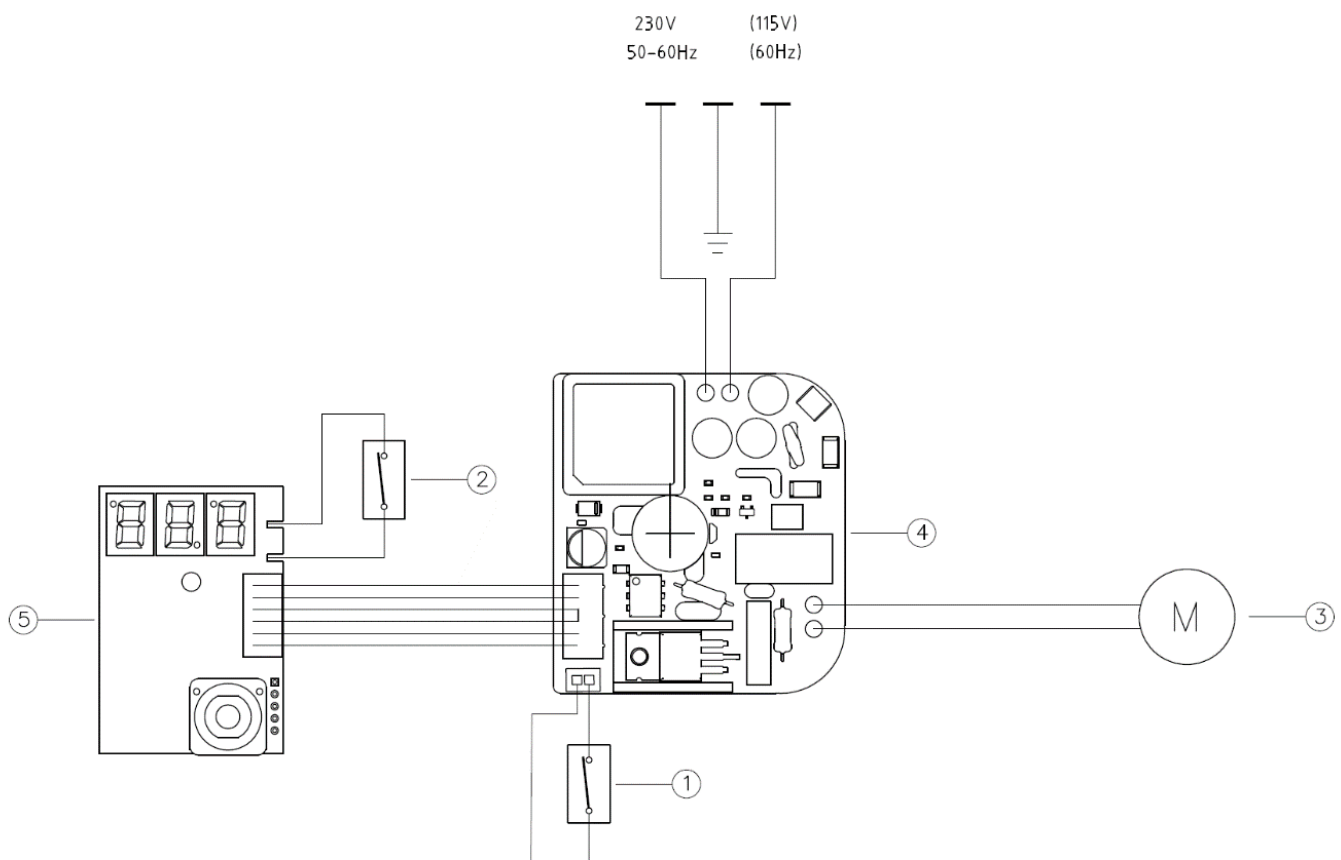
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of the following directives:

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2006/42/EC Machinery directive

2014/30/EU Electromagnetic compatibility directive

2015/863/EU (RoHS III) Restriction of the use of certain hazardous substances

2012/19/EU (WEEE) Waste of electric and electronic equipment

plus modifications / più modifiche / plus modifications / más sucesivas modificaciones / in der jeweils gültigen Fassung /
加修改.

12. Declaration of conformity ^{UK}_{CA}

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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VELP Scientifica Srl