




MEDIA INLET LINE

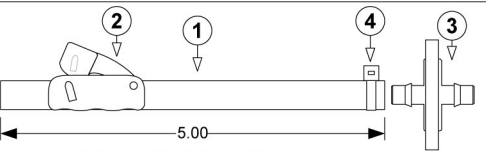


SPARGER LINE

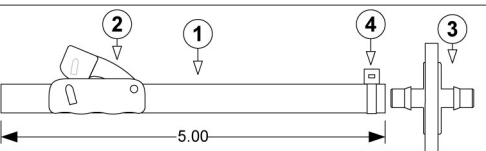


OVERLAY LINE

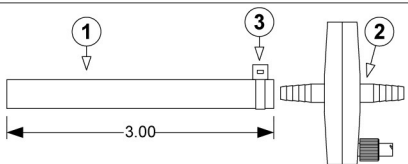

CONDENSER / VENT LINE

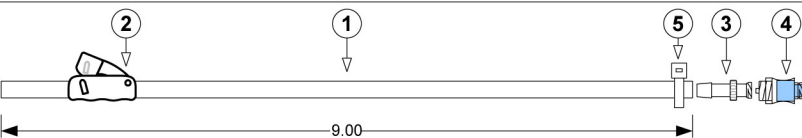

SAMPLE LINE

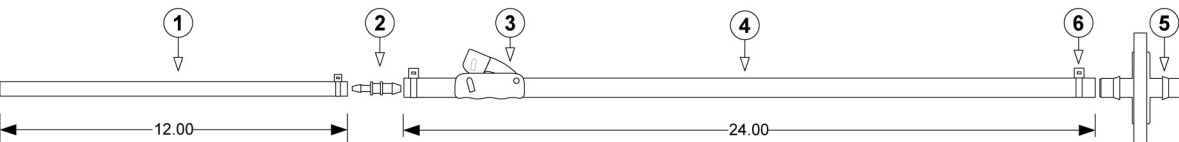

TRI-PORT LINE

Material List

ITEM	QTY	DESCRIPTION
MEDIA INLET LINE		
1	1	.1875" ID x .375" OD PLATINUM CURED SILICONE
2	1	PURE FIT TUBE CLAMP, POLYPRO
3	1	25 mm 0.2µm GAMMA STABLE HYDROPHOBIC FILTER
4	1	TIE WRAP, NYLON
SPARGER LINE		
1	1	.1875" ID x .375" OD PLATINUM CURED SILICONE
2	1	PURE FIT TUBE CLAMP
3	1	25 mm 0.2µm GAMMA STABLE HYDROPHOBIC FILTER
4	1	TIE WRAP, NYLON
OVERLAY LINE		
1	1	.1875" ID x .375" OD PLATINUM CURED SILICONE
2	1	PURE FIT TUBE CLAMP
3	1	25 mm 0.2µm GAMMA STABLE HYDROPHOBIC FILTER
4	1	TIE WRAP, NYLON
CONDENSER / VENT LINE		
1	1	.1875" ID x .375" OD PLATINUM CURED SILICONE
2	1	50 mm 0.2µm GAMMA STABLE HYDROPHOBIC FILTER
3	1	TIE WRAP, NYLON
SAMPLE LINE		
1	1	.063" ID x .188" OD 374 C-FLEX
2	1	PURE FIT TUBE CLAMP, POLYPRO
3	1	FEMALE LUER x .0937" HOSE BARB, POLYPRO
4	1	SWABABLE LUER VALVE, POLYCARBONATE
5	1	TIE WRAP, NYLON
TRI-PORT LIQUID LINE		
1	1	.0625" ID x .1875" OD PLATINUM CURED SILICONE
2	1	.125" x .0937" REDUCING CONNECTOR, POLYPRO
3	1	PURE FIT TUBE CLAMP, POLYPRO
4	1	.125" ID x .250" OD 374 C-FLEX
5	1	25 mm 0.2µm GAMMA STABLE HYDROPHOBIC FILTER
6	3	TIE WRAP, NYLON